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Using JT 65T adjustable inserts in conjunction with optic holders 65,  $\varnothing 45$  mm or cubes W 65, highly stable beam deflectors can be assembled. With cubes 65, it is a simple matter to install lamps up to 50 W and condenser mirrors, or build a microscope set-up with two viewing points. By combining several cubes, precise ring-resonators, and Mach-Zehnder or Michelson interferometers can be built.

The adjustable elements can be fitted to any side of a cube using M3 screws, or to optic holders 65 with  $\varnothing 45$  mm fixing hole.

In the adjustable inserts T the optics holder is supported on packets of disc springs against the flange component, and can be adjusted with 3 screws. They are very precise, shock and vibration resistant, and stable over long periods. The third screw provides a limited possibility to compensate for different substrate thicknesses, or to eliminate reflections.

The adjustable mount 65T for plane mirrors is designed so that it can be installed in a cube 65 together with a second adjustable mount, for example a beam splitter.

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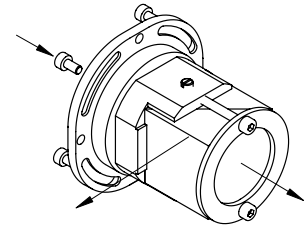
### Adjustable Insert JT 65T for Beam Splitter Cube and Prisms

for elements with 30 mm edge length. Lateral mounting means that prisms can also be installed. Used in interferometry, it is suitable for beam diameters up to 25 mm.


		Typ/type	Bestell-Nr./part no.
Justiereinsatz JT 65T für Strahlteilerwürfel oder Prisma 30x30x30 mm	adjustable insert JT 65T for beam splitter cube or prism 30x30x30 mm	JT 65T-STW30Z-O	16.214.1030
Justiereinsatz JT 65T mit Strahlteilerwürfel STWU-30-BK7-L/4-50:50, Bestell-Nr. 60.421.3050, 30x30x30 mm	adjustable insert JT 65T with beam splitter cube STWU-30-BK7-L/4-50:50, part no. 60.421.3050, 30x30x30 mm	JT 65T-STW30Z-M	16.214.1130

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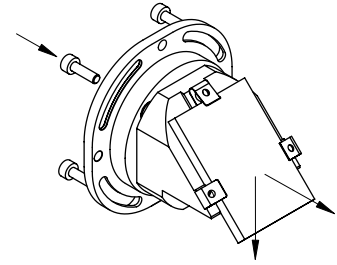### Adjustable Insert JT 65T for Plates

for adjusting plates, e.g. beam splitter plates, 30x42.5x4 mm thick. The plates are retained by springs at three points. The stress-free mounting means that the insert can be used for interferometry, even at large beam diameters.

Justiereinsatz JT 65T für Strahlteilerplatte 30x42,5x4 mm	adjustable insert JT 65T for beam splitter plate 30x42.5x4 mm	JT 65T-STPZ-O	16.214.2030
Justiereinsatz JT 65T mit Strahlteilerplatte STPL-30x42.5-BK7-L/4-50:50, Bestell-Nr. 60.411.3050, 30x42,5x4 mm	adjustable insert JT 65T with beam splitter plate STPL-30x42.5-BK7-L/4-50:50, part no. 60.411.3050, 30x42.5x4 mm	JT 65T-STPZ-M	16.214.2130

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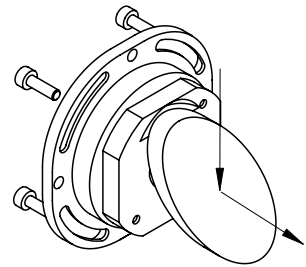
Optimale Spiegeldicke: 4 mm.

### Adjustable Insert JT 65T for Deflection Mirror

for deflecting expanded beams at 90°. Stress-free bonding to the optical glass substrate makes this suitable for high-quality laser mirrors. For beam diameter up to 30 mm suitable.

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
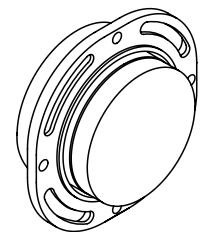
### Adjustable Insert JT 65T for Mirrors with Flat Rear Surface

Stress-free bonding to an optical glass substrate makes this suitable for reference mirrors in interferometers. The biggest diameter of the mirror with complete flexibility is 40 mm.

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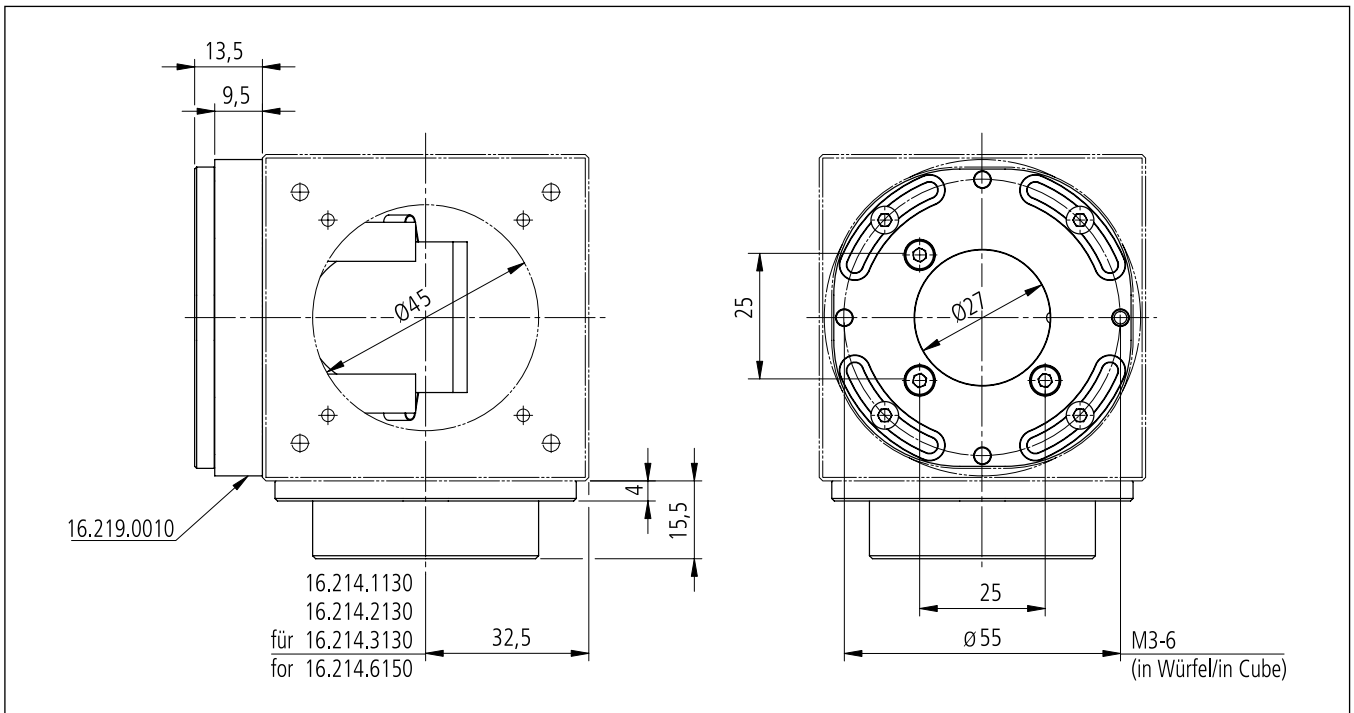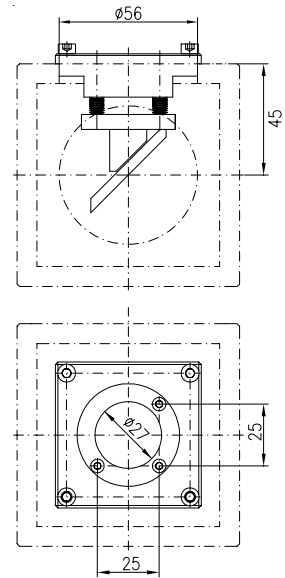
## Adjustable Insert for Mounting Cube 65-4

These adjustable inserts are specially adapted for installation in mounting cubes 65-4. In all other respects they are identical to adjustable inserts JT 65.

incl. 4 screws ISO 4762 M4x8 and  
4 threaded inserts M8

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